

Human Topoisomerase I Microplate Assay Kit



Product Description (Product Numbers TRT101 and TRT102)

The kit is supplied with sufficient human topo I enzyme, plasmid DNA substrate (supercoiled pNO1; supplied at 1 mg/ml), 10X Assay Buffer, Enzyme Dilution Buffer and TFO1 oligo for 100 assays. The enzyme is supplied at a concentration of 10 U/μl in Dilution Buffer. The kit is also supplied with sufficient Wash Buffer, TF buffer and T10 buffer for one 96-well plate. These buffers are supplied as 20X concentrates and must be diluted prior to use with ultra-pure water.

Store enzyme at -80 °C. (Stable for 3 months undiluted.) It is recommended that the enzyme is aliquoted to avoid repeated freeze-thaw cycles.

For *in vitro* laboratory research use only.

Dilution Buffer

10 mM Tris.HCl (pH 7.5)
1 mM DTT
1 mM EDTA
50 % (v/v) glycerol
100 μg/ml albumin

Assay Buffer (supplied as 10X stock)

20 mM Tris.HCl (pH7.5)
200 mM NaCl
0.25 mM EDTA
5 % (v/v) glycerol

TF Buffer (supplied as a 20X stock)

50 mM sodium acetate (pH 5.0)
50 mM NaCl
50 mM MgCl₂

Wash Buffer (supplied as a 20X stock)

20 mM Tris.HCl (pH 7.6)
137 mM NaCl
0.01 % (w/v) BSA
0.05 % (v/v) Tween-20

T10 Buffer (supplied as a 20X stock)

10 mM Tris-HCl (pH 8)
1 mM EDTA

Preparation of Plate and Relaxation Assay

Rehydrate wells with 3 x 200 μl Wash Buffer (diluted from 20X stock before use).


Immobilize 100 μl of 500nM TFO1 oligo in well (5 μl of 10 μM TFO1 in 95 μl Wash Buffer), 5 minutes at room temperature. Wash off excess oligo with 3 x 200 μl Wash Buffer.

Incubate 1.5 U of human topo I with 0.75 μg of supercoiled pNO1 in a reaction volume of 30 μl at 37°C for 30 minutes in Assay Buffer. Incubate reaction in well.

Add 100 μl TF Buffer (diluted from 20X stock before use) to well and incubate for a further 30 minutes at room temperature to allow triplex formation.

Remove liquid from well and wash with 3 x 200 μl TF Buffer to remove unbound plasmid.

Stain with appropriate fluorescence stain (Suggested stain, SYBR Gold® (Invitrogen) diluted to 1X with T10 buffer. Add 200 μl per well. Incubate for 10 - 20 minutes, mix and read in fluorescence plate reader; Ex: 495 nm; Em: 537 nM).

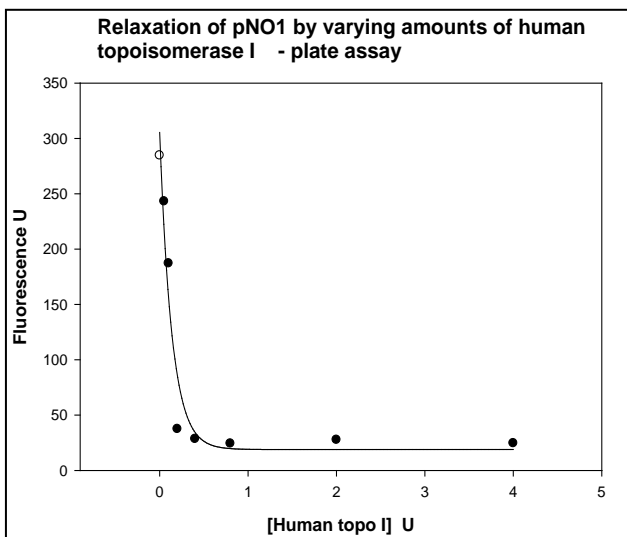

Quality Control

1) Purity: Human topoisomerase I is purified to > 95 % purity as judged by SDS-polyacrylamide gel electrophoresis. 2) Tests for human topoisomerase II contamination by looking for decatenation of kDNA under topoisomerase II assay conditions were negative. 3) kDNA or pBR322 were also incubated for 4hrs in assay buffer + 10 mM MgCl₂ at 37 °C. These tests were negative for the formation of linear products, showing the absence of nuclease contamination.

References

Maxwell, A., Burton, N.P. and O'Hagan, N. (2006) High-throughput assays for DNA gyrase and other topoisomerases. *Nucleic Acid Res.* **34(15)**, e104

Patent held by Plant Biosciences Ltd., Norwich, Norfolk, UK (Patent No. GB0424953.8.)

Kit issued with limited licence for individual use only